



Kentek Laser Safety Containment System Curtains and Barriers

Use Maintenance, and Cleaning

General Use

The primary use of your Laser Safety Containment System is to provide a safe environment in which to operate the laser. Follow installation instructions carefully and contact your Kentek representative for assistance.

Maintenance

Perform a routine inspection of the Laser Safety Containment System to verify that it functions as intended. Inspect curtains and barriers for tears or holes in the fabric and any other potential hazard. If you find an area of concern, discontinue use and contact your Kentek representative. Your Kentek representative can suggest options for repair of the area or other solutions for the problem.

Cleaning

Kentek's FLEX-GUARD® laser safety curtains can be cleaned with the following cleaning agents:

- A mild detergent solution, such as Dawn.
- 30% bleach
- 100% isopropyl alcohol
- Super Sani-Cloth® Germicidal Disposable wipes
- Less than 3% peroxide solution

To prevent damage to the laser safety curtain, use a soft bristle brush.